

Quick Reference Guide for Vented Gas Cabinet for Cylinders

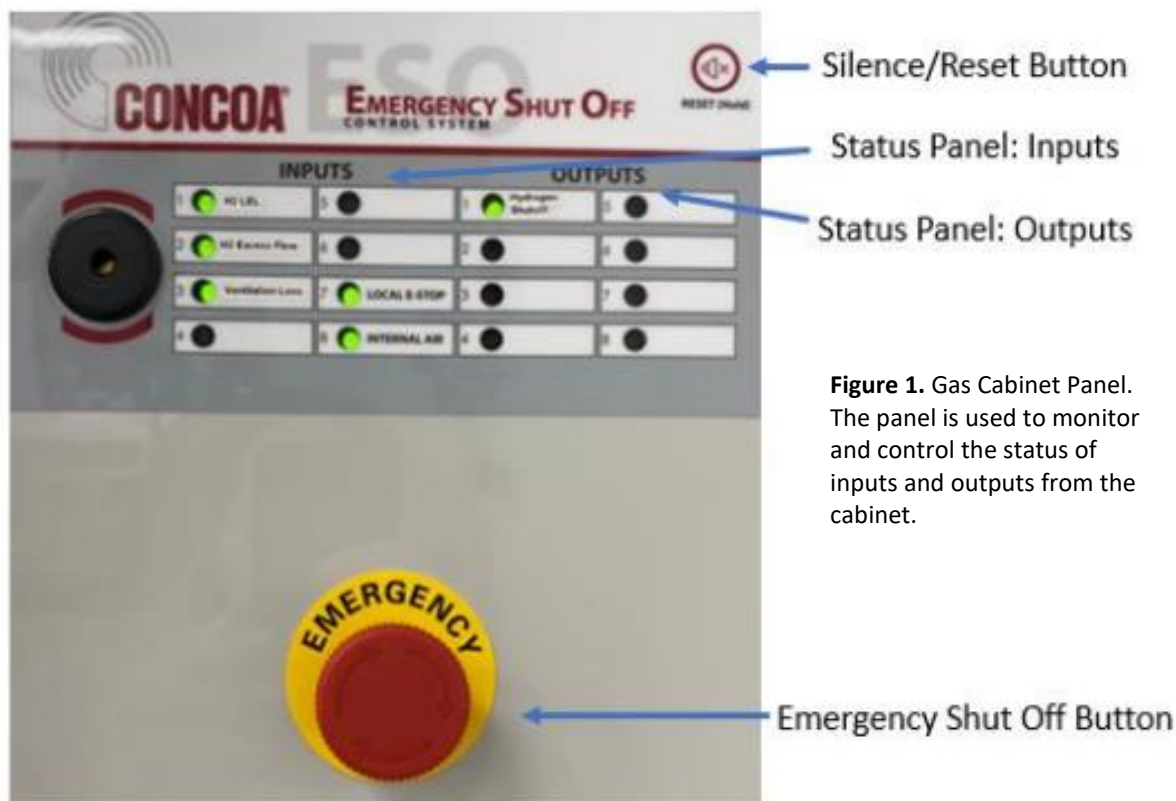


Figure 1. Gas Cabinet Panel. The panel is used to monitor and control the status of inputs and outputs from the cabinet.

In Case of Emergency:

In case of emergency, press the Emergency Shut Off Button. This will shut off the gas. Afterwards, please contact EHS at 410-704-2949 or safety@towson.edu or the Building Manager (contact listed below) to alert them to the situation.

How to shut off alarm:

If alarm is sounding, push the Silence/Reset Button in the top right corner. This will not turn the gas back on, but it will silence the noise.

How to reset:

If the Status Panel has all lights green (steady or flashing) on the Input column, push and hold the Silence/Reset Button for 5-10 seconds until the Outputs column light turns green; the gas is now able to be used. If the Inputs column has a red light, please contact the Building Manager for assistance.

SC Building Manager: Josh Wilhide, 410-704-5229